

# xxsealer™

PCR plate heat sealer with programmable temperature and dwell



## User Manual



**BJS BIOTECHNOLOGIES**  
A LIFESCIENCE COMPANY

**This Manual Copyright © 2014 BJS Biotechnologies**

All rights reserved. No part of this manual may be reproduced, stored in a retrieval system, or transmitted, by any means or in any form, without the prior permission of BJS Biotechnologies Ltd (BJS Bio).

The information contained in this manual is subject to change without notice. BJS Bio assumes no responsibility for any errors that may appear in this or related documentation.

Document: XP-5003-01-1  
Revision: 1.0 Jan 2014

**xxpress® xxplates, and xxsealer** are the brand names for BJS Biotechnologies own range of PCR thermal cycler products and accessories. For further details please visit our web site.

[www.xxpressPCR.com](http://www.xxpressPCR.com)

All sales and technical enquiries should be addressed to:

BJS Biotechnologies Ltd  
65 Bideford Avenue  
Perivale  
Greenford  
Middlesex  
UB6 7PP  
United Kingdom

+44 (0)203 021 3750

[sales@xxpressPCR.com](mailto:sales@xxpressPCR.com) or [support@xxpressPCR.com](mailto:support@xxpressPCR.com)

## Contents

1	Glossary of Terms and Abbreviations .....	4
2	Symbols Used in this Instruction Manual .....	5
3	Unit Description .....	6
4	Technical Specification.....	9
5	Safety Precautions and Limitations of Use.....	10
6	Unit Installation .....	14
7	Unit Operation .....	16
7.1	User Display and Controls.....	16
7.2	Operating the Controls of the xxsealer.....	17
7.3	Performing an xxplate Sealing Operation.....	18
7.4	Switching Off the Audible Alert.....	20
8	Maintenance and Servicing .....	21
8.1	Replacing the Unit Fuse .....	21
8.2	Over Temperature Safety Cut-out .....	22
8.3	Routine Cleaning and Inspection .....	22
8.4	Decontamination Procedure .....	23
8.5	Transportation and Storage.....	23
9	Warranty and Returns .....	25
10	Product Disposal .....	26

## Tables









<b>Table 1: Advisory Symbol Meanings.....</b>	<b>5</b>
<b>Table 2: Unit Features.....</b>	<b>6</b>
<b>Table 3: Plate and Adaptor Block Types.....</b>	<b>7</b>
<b>Table 4: Recommended Film and Foil Types .....</b>	<b>8</b>
<b>Table 5: Mains Voltage Selection.....</b>	<b>14</b>
<b>Table 6: User Buttons and Status LEDs.....</b>	<b>16</b>

## 1 Glossary of Terms and Abbreviations

<b>ANSI</b>	American National Standards Institute
<b>BJS Bio</b>	BJS Biotechnologies Limited, UK Company Number 02959160
<b>Deepwell Plate</b>	Plate with an SBS footprint featuring 48, 96 or 384 wells with a larger volume than microplates
<b>DWP</b>	Deepwell plate
<b>EMC</b>	Electro-Magnetic Compatibility
<b>Microtitre Plate</b>	Plate with an SBS footprint featuring 24, 48, 96 or 384 wells
<b>MTP</b>	Microtitre plate
<b>PCR</b>	Polymerase Chain Reaction
<b>PPE</b>	Personal Protective Equipment
<b>SBS</b>	Society for Bio molecular Screening
<b>Semi-skirted PCR Plate</b>	PCR plate with an outer surrounding half edge
<b>Skirted PCR Plate</b>	PCR plate with an outer surrounding edge
<b>Un-skirted PCR Plate</b>	PCR plate without an outer surrounding edge
<b>Well</b>	A single sample cavity in a Microtitre plate, PCR plate or Deepwell plate
<b>xxplate</b>	BJS Bio PCR test plate for the xxpress real-time thermal cycler. Available in 24, 54, and 96 well formats.

## 2 Symbols Used in this Instruction Manual


The following advisory symbols are used in this manual.

Table 1: Advisory Symbol Meanings	
 <b>DANGER</b>	Indicates a Risk of Electric Shock which could, if not avoided, result in severe injury or death.
 <b>DANGER</b>	Indicates a Burn Hazard which could, if not avoided, result in severe injury or death.
 <b>DANGER</b>	Indicates a Risk of Explosion which could, if not avoided, result in severe injury or death.
 <b>WARNING</b>	Indicates a hazardous situation which could, if not avoided, result in severe injury or death; or severely damage the unit.
 <b>CAUTION</b>	Indicates a hazardous situation which could, if not avoided, result in minor or moderate injury; or degrade or impair the functionality of the unit.
 <b>CAUTION</b>	Indicates a Risk of Crush hazard due to moving parts which could, if not avoided, result in minor or moderate injury.
	Advisory note or other useful information.
	Refer to "Section NN" for more details.

### 3 Unit Description

The **xxsealer** provides a safe and controlled method for sealing plastic sample plates to protect the samples from evaporation and contamination during the heating and cooling of a PCR process. The sealer is adjustable for both the sealing time and temperature, and gives a visual countdown display during the sealing process for consistent results.





The unit has the following external features:

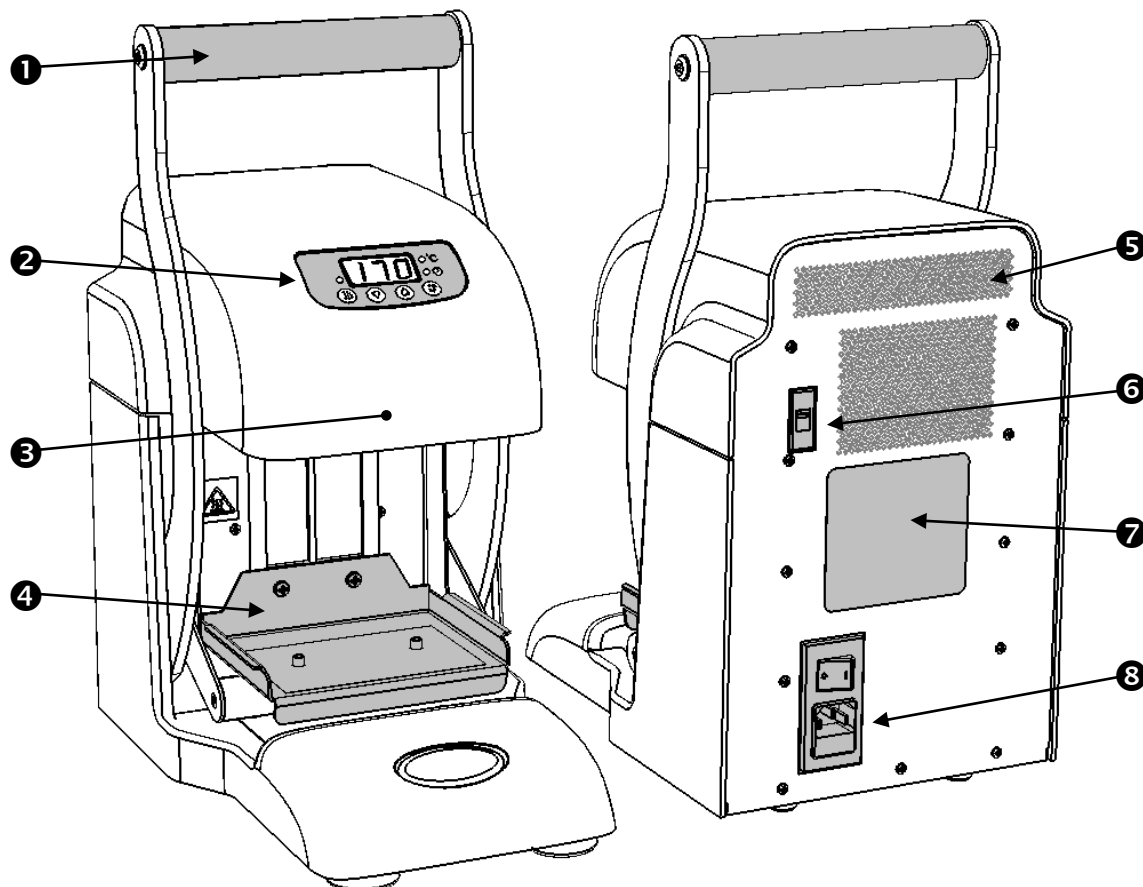

**Table 2: Unit Features**

①	Seal Activation Handle	⇒ 7.2
②	User Interface Display and Control Buttons	⇒ 7.1
③	Heater Plate (Internal)	⇒ 7.1
④	Adaptor Block Carrier	⇒ <b>Table 3</b>
⑤	Rear Ventilation Holes	⇒ 5
⑥	Mains Voltage Selector	⇒ 5
⑦	Product Information Label	⇒ 4
⑧	Power Inlet, On/Off Switch and Fuse Holder	⇒ 5

The **xxsealer** unit is designed to be used with a range of sample plates and consumables – some of which are listed below:
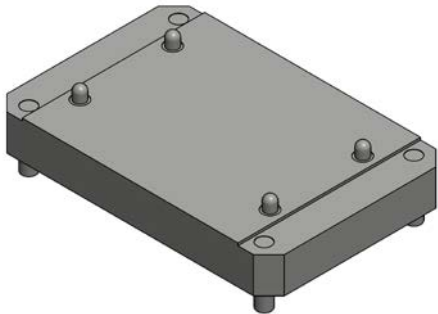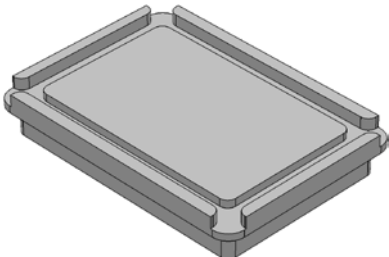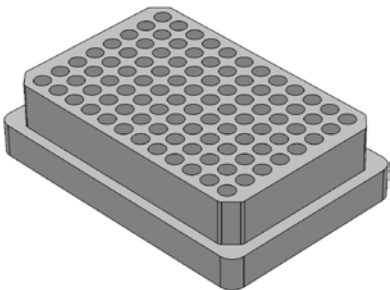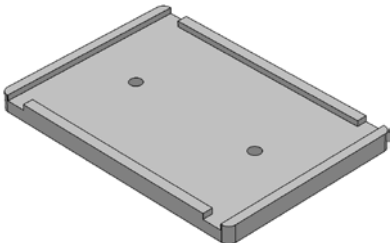
<b>Table 3: Plate and Adaptor Block Types</b>	
<b>Adaptor Block Type</b>	<b>Suitable Plate Types</b>
	<b>BJS xxplate</b> (Suitable for 24, 54, and 96 well xxplates)
	96 Well Microtiter Plate 384 Well Microtiter Plate PCR 96 Well Skirted Plate
	PCR 96 Well Semi-skirted Plate PCR 96 Well Un-skirted Plate
	Deepwell Plates
<b>Plate Material Types</b>	Polypropylene, Polyethylene or Polystyrene

Other specialist adaptor plates may be available at request. Please contact your distributor for details.

**Table 4: Recommended Film and Foil Types**

<b>Table 4: Recommended Film and Foil Types</b>	
<b>Films</b>	Clear Polyester/Polypropylene Laminate Clear Polymer Thin Clear Polymer
<b>Foils</b>	Foil Polypropylene Laminate Foil Laminate Foil

**Note:**

The only film approved for use with **xxplates** is the BJS Bio's **xxplate** sealing film. Part number HS-3001-01-1.





## 4 Technical Specification

### Model Type

Model Name	<b>xxsealer</b>
Model No	VTS-953-A

### Physical Unit Properties

Dimensions (W x D x H)	220 mm x 325 mm x 425 mm
Weight (without adaptor)	7.2 kg

### Mains Supply

Power Cord Rating	IEC C13, 3-Core, 5A min
Inlet Module Type	IEC C14, DPST, Single Fuse
Supply Voltage Selections	<input type="checkbox"/> 115 110 to 120 VAC $\pm 10\%$ <input type="checkbox"/> 230 220 to 230 VAC $\pm 10\%$
Supply Frequency Range	50 to 60 Hz $\pm 5\%$
Power Consumption	350 W max
Fuse Ratings and Sizes	<input type="checkbox"/> 115 T3.15AH 250V 20x5mm <input type="checkbox"/> 230 T1.6AH 250V 20x5mm

### Operating Environment

Temperature Range	+18 to +30 °C
Relative Humidity Range	20% to 80% non-condensing
Maximum Operating Altitude	2200 m above sea-level

### Storage and Transportation

Temperature Range	-10 to +50 °C
Relative Humidity Range	20% to 95% non-condensing

### Sealing Parameters

Temperature Range	125 to 200 °C in 1 °C steps
Timer Range	1.0 to 9.0 seconds in 0.5 second steps
Audible Alert	On or Off

## 5 Safety Precautions and Limitations of Use

It is essential that all users of this equipment have fully read and understood the following safety precautions and limitations of use before installing or operating the **xxsealer** unit.

### IMPORTANT



**WARNING**

The protection provided by this equipment may be impaired if it is not used in a manner described in this manual.



**WARNING**

It is essential that the user of this equipment is aware of the potential hazards associated with the unit and its accessories. All operators should be familiar with the safety precautions and warnings given in these instructions before attempting to operate the unit. Improper use of this unit or its accessories may impair their functionality and invalidate the manufacturer's warranty.

### Unit Handling Precautions



**CAUTION**

Care should be taken not to drop the unit or subject it to rough physical handling, both during normal use and during transportation and storage.



**WARNING**

The unit should be held by the handle when being lifted or moved. Do not lift the unit by any other part of the casework. Care should be taken when lifting the unit due to its weight.



**CAUTION**

Care should be taken not to knock the LED display. Do not use excessive force when pressing the control buttons or operating the seal activation handle.

## Unit Installation and Operating Environment



**DANGER**

The **xxsealer** unit is designed for indoor laboratory use only. The acceptable operating temperature range is 18°C to 30°C, with a relative humidity of 20% to 80% non-condensing, at a maximum altitude of 2200m above sea level.



**WARNING**

If the unit is stored in conditions outside of these ranges, it must be left to stand unpowered until it has acclimatised to within these environmental limits before being powered.



**DANGER**

Use only the AC mains power cord provided with the unit or as specified Section 4.

The unit must be connected to a suitably earthed mains supply, with appropriate earth-leakage and over-current protection.



**WARNING**

Always ensure that the mains power connector is securely inserted into the rear of the unit, and any excess power cord does not pose a potential trip or pull hazard.



**DANGER**

Do not operate the unit in any area which is, or has been, or is thought to have been exposed to explosive or flammable gases, vapours or liquids.



**WARNING**

The unit must be installed and operated on a solid, stable and level working surface; ensuring that the ventilation holes at the rear of the unit are not obstructed.

## General Operating Precautions



**DANGER**

Ensure that the power is switched off at both the AC mains supply outlet and at the back of the unit before inserting or removing the mains power cord.



**DANGER**

The heater plate can reach temperatures of 200°C and will remain hot for a considerable time after being turned off.

Extreme care must be taken not to touch the heater plate as it will cause a severe burn injury.



**DANGER**

The unit is intended for use with plates containing biological samples only.

Never use the unit to seal any explosive, volatile or highly reactive substances or chemicals.



**CAUTION**

There is a possible finger crush hazard due to the moving parts of the handle and plate carrier.

Care should be taken when operating the handle.

## Unit Maintenance and Serviceability



**DANGER**

There are no user or operator serviceable parts inside the unit.  
Do not remove the unit casework.



**WARNING**

Removal of the unit's casework will void the manufacturer's warranty and may expose the user to a Risk of Electric Shock resulting in serious injury or death.



**DANGER**

Once installed, the externally accessible unit fuse will only blow under a fault condition. This fuse should only be changed after the unit has been thoroughly inspected by a qualified engineer. See Section 8.1 for details.



**DANGER**

Always switch off the unit and disconnect the power cord before performing any cleaning or decontamination procedure.  
If liquid is spilt into or over the unit, switch off and disconnect the power from the AC mains outlet before attempting to deal with the spillage.



**CAUTION**

Ensure that the heater plate has cooled down to room temperature before performing any cleaning operation and before moving or storing the unit.



**CAUTION**

The use of harsh chemicals and cleaning agents may damage the unit and degrade its performance.  
Always follow the cleaning and decontamination procedures specified in Sections 8.3 and 8.4 of this instruction manual.



**CAUTION**

Do not autoclave any part of the unit or its accessories.

## 6 Unit Installation

Before installing the **xxsealer** unit, please check that the delivery is complete and that the unit and any accessory parts are intact and free from any signs of transportation damage. Also ensure that all external and internal packaging has been removed from the unit before installation.



Please retain all packaging for future transportation and storage of the unit and its accessories.

The **xxsealer** unit should be installed in a location which meets the following requirements:



- Safe and suitable operating environment (see Section 5)
- Solid, stable, level working surface
- At least 10cm clearance around the unit to adjacent objects and walls
- Earthed AC mains power connection (see Section 4)



**WARNING**

Please also observe and abide by the **Unit Installation and Operating Environment** safety precautions and preconditions listed in Section 5.



It is important to check that the Mains Voltage Selector switch on the rear of the unit has been set to the correct position for the desired operating voltage, and the correct value of fuse has been fitted (see Table 5).

<b>Table 5: Mains Voltage Selection</b>		
<b>Operating Voltage</b>	<b>110 to 120 VAC</b>	<b>220 to 230 VAC</b>
<b>Voltage Selection</b>	115 	230 
See Section 3 for switch location (Ⓔ) on unit rear panel.		
<b>Fuse Rating</b>	3.15AH	1.6AH



**WARNING**

Ensure that the correct fuse type has been fitted in the mains inlet fuse holder for the desired voltage selector switch value. See Table 5 for details.

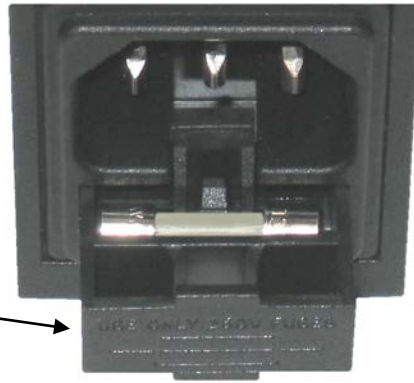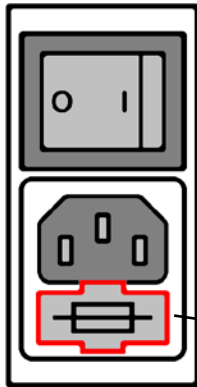

1. Remove the AC plug from the power inlet module
2. Use a small flat bladed screwdriver to carefully pull out the fuse holder
3. Fit the correct fuse in the far position only
4. Push the holder back into the inlet mode

Install and test the **xxsealer** unit using the following procedure:

- 1) Place the unit on the suitably selected working surface (as specified above), ensuring that the ventilation holes on the rear of the unit are not obstructed.
- 2) Connect the unit to the AC mains power outlet using the mains power cord supplied or as specified in Section 4.
- 3) Switch the mains power on at supply outlet first, and then switch the unit on using the power switch located at the rear of the unit.
- 4) Check that the unit is stable and safe by performing a trial sealing operation, as described in Section 7.2.



**DANGER**

If the unit has been stored in a cool environment, it must be left to stand unplugged until it has acclimatised to the new room temperature before being powered.

## 7 Unit Operation




**WARNING**







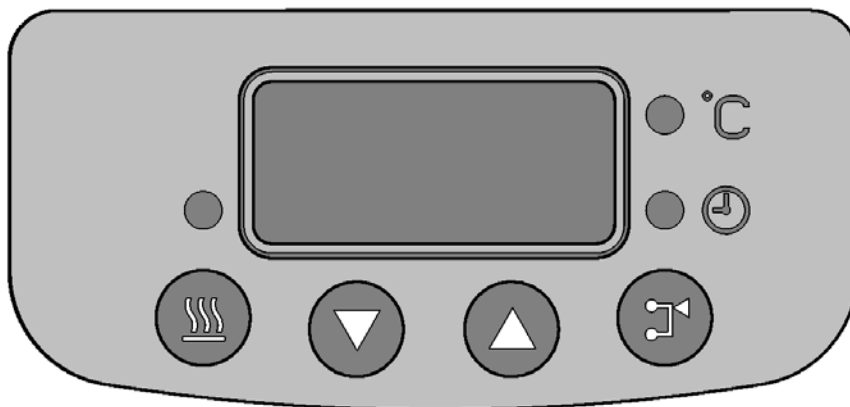



Please ensure that you have read and fully understood all of the **Safety Precautions and Limitations of Use** listed in Section 5 before attempting to operate the **xxsealer** unit.

### 7.1 User Display and Controls

The unit's user interface consists of a large 3-digit LED display, four control buttons and three status indicator LEDs.



The LED display shows the current temperature or sealing time set-point (depending on the mode). The button and LED functions are listed in Table 6.

Table 6: User Buttons and Status LEDs		
 	<b>Heater Status LED</b> <b>HEATER On/Off Button</b>	 <b>Off</b> – Heater Turned Off  <b>Flashing</b> – Warming Up  <b>On</b> – Unit Ready For Use
	<b>DOWN</b> Button	
	<b>UP</b> Button	
	<b>MODE</b> Select Button	
	<b>Temperature</b> Mode LED	
	<b>Timer</b> Mode LED	



## 7.2 Operating the Controls of the xxsealer

When the unit is switched on, it recalls the last sealing operation settings and the displays the sealing temperature or time (depending on the display mode), and automatically restores the heater on/off state.

Pressing the **MODE** button toggles the LED display between the plate heater **Temperature** set-point in degrees centigrade and the seal **Timer** setting in seconds.

The sealing temperature is set by pressing the **MODE** button to illuminate the **Temperature** LED (● °C) then using the **UP** and **DOWN** buttons to set the temperature in the range 125 to 200 °C in 1°C steps.

Likewise, the sealing time is set by pressing the **MODE** button to illuminate the **Timer** LED (● ⌚) then using the **UP** and **DOWN** buttons to set the time in the range 1 to 9 seconds in 0.5 second steps.

Pressing the **HEATER** button toggles the plate heater between on and off. When the **Heater** LED is flashing, the heater is on and the plate is warming up, which will normally take about 10 minutes.

Once the **Heater** is on the LED display shows the current temperature or sealing time set-point (depending on the mode).

When the heater plate is up to temperature (within 2°C of the set-point) the **Heater** LED will remain constantly on, and the **xxsealer** is ready to begin sealing.

The operator may set a precise temperature and seal time for the sealing operation. Sealing is initiated by pulling the handle down, which raises the sample plate into contact with the heater plate. The unit detects when the sample plate is correctly positioned and starts a countdown timer. When the display reads zero, the operator moves the handle back to its upmost position, which moves the sample plate away from the heater plate. The sample plate is then manually removed for storage or further processing.


### 7.3 Performing an xxplate Sealing Operation

A typical operating sequence to seal an **xxplate** is:

- 1) Switch the unit on and set the sealing time to 1.5 seconds and the heater plate set-point temperature to 170°C (using the method described above). These are the recommended setting for the **xxplate** film.
- 2) If necessary, press the **HEATER** button to turn the heater on.
- 3) Allow sufficient time for the unit to reach temperature (the **Heater** LED will stop flashing and remain constantly on).
- 4) Fit the **BJS xxplate Adapter Block** into the **PCR 96 Well Skirted Plate Adapter Block** (both provided with the **xxsealer**) and locate in the **Adapter Block Carrier**




- 5) Load the **xxplate** onto the **xxplate Adapter Block** locating it onto the sprung pins. It should sit flat on the adapter block. And add the sealing film on top, the correct way up.

	Care should be taken not to touch the surface of the heater plate whilst loading the sample plate.
WARNING	


6) Grasp the handle with one or two hands, as appropriate, and pull the handle forward and down to raise the plate carrier until the sample plate contacts the heater plate.



7) When correct sealing pressure is achieved, an audible alert will sound (if enabled – see Section 7.4) and the timer will start counting down from the set value. When zero seconds is reached, the display will flash and the audible alert sound.

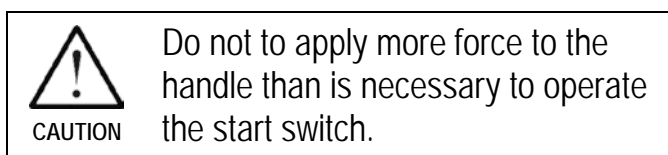

8) Immediately return the handle to its upmost position to retract the sealed sample plate from the heater plate.

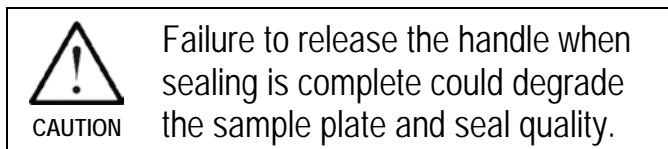

9) Remove the sealed plate from the unit and review the seal integrity.



For optimum performance, the sealing temperature and/or time may need to be adjusted to best suit the sample plate and film/foil types being sealed. Various types of plastic sample plate may be loaded onto a plate carrier and thermal sealed using a suitable foil or clear film. See Section 3 for details.

## 7.4 Switching Off the Audible Alert



The unit is fitted with an audible alert which sounds at the beginning and end of each sealing operation.

If desired, this alert can be turned off (or back on again) using the following procedure:

- 1) If necessary, press the **MODE** button to select the **Temperature** LED setting mode.
- 2) Next, simultaneously press both the **HEATER** and **MODE** buttons together.
- 3) The display should show “01” or “00”. If it does not, repeat this procedure from step 1 above.
- 4) Now press the **DOWN** button to turn the alert off (display shows “00”) or the **UP** button to turn the alert back on (display shows “01”).
- 5) Finally, press the **MODE** button to store the new setting.

## 8 Maintenance and Servicing


Although the **xxsealer** unit does not require any scheduled maintenance or servicing, the operator should regularly clean and inspect the unit for any detects, as described in Section 8.3 below.

	<p>Please observe and comply with all of the <b>Unit Maintenance and Serviceability</b> precautions listed in Section 8.</p>
<p><b>DANGER</b></p>	<p>Never removed the unit casework. There are no user or operator serviceable parts inside the unit.</p>
	<p>Always switch off and unplug the unit before performing any cleaning or disinfecting tasks.</p>
<p><b>WARNING</b></p>	


For technical and service related enquiries, please contact your distributor or BJS Bio at the address given on Page 2 of this manual.

### 8.1 Replacing the Unit Fuse

The unit fuse should only be replaced by a suitably qualified technician.

	<p>The unit fuse will only blow as a result of an internal unit fault or if the voltage selector switch has been incorrectly set (see Section 5). This fuse should only be changed after the unit has been thoroughly inspected, and must be replaced with the exact type specified in Section 4.</p>
<p><b>DANGER</b></p>	

Thoroughly inspect the unit for any signs of damage, loose components or liquid spillage or ingress. If in doubt, please contact BJS Bio on the number given on Page 2 of this manual.

	<p>The fuse holder is removed by disconnecting the mains cord and then using a small flat bladed screwdriver to carefully pull open the fuse holder (see Section 5) and remove the old fuse.</p> <p>After replacing the fuse with the corrected rated one for the operating voltage being used (see Table 5), push the fuse holder firmly back into the inlet module.</p> <p>The unit must be electrically safety tested for excess leakage current before being repowered from the mains supply.</p>
<p><b>DANGER</b></p>	

## 8.2 Over Temperature Safety Cut-out

The unit is fitted with a non-resettable thermal fuse which blows if the heater plate temperature exceeds 235 °C.

In the unlikely event of a fault condition, this fuse will permanently disable the heater to protect the user and unit from injury and damage. If this occurs, the heater plate will remain cold and the **xxsealer** will need to be returned for repair.

## 8.3 Routine Cleaning and Inspection

The unit casework should be cleaned and inspected at regulator internals, and whenever contamination or spillage occurs, as follows:

1. Switch off the unit and disconnect the power before performing any inspection checks or cleaning.
2. Before cleaning, always inspect the unit casework, heater plate and moving parts for any signs of wear, damage, cracks or other defects.
3. Clean the casework using a damp cloth soaked with a disinfectant solution (such as Virkon), whilst wearing suitable PPE.
4. Carefully clean the heater plate surface to remove any debris or sealing material.
5. Remove any debris or fluff from around or between the moving parts of the handle mechanism and sliding plate carrier.
6. Clean the display and buttons, taking care to avoid over wetting.
7. Check that the ventilation holes on the rear of the unit are clear of dust and fluff build-up.
8. Check and clean the adaptor blocks.



**DANGER**

After cleaning, ensure that the unit is thoroughly dry, especially around the mains power inlet, before reconnecting the power cord and switching the unit on.

## 8.4 Decontamination Procedure

The unit and accessories should be decontaminated using the following procedure before being stored or transported.

### Decontamination Procedure

Thoroughly clean all outside surfaces of the product (including any accessories, power cords, manuals, packaging, etc) with a damp cloth soaked with suitable disinfectant solution (such as Virkon).

## 8.5 Transportation and Storage

The **xxsealer** unit and its accessories should be thoroughly decontaminated using the procedure detailed in Section 8.4 before being placed in its original packaging for transportation or storage.



### WARNING

Refer to Section 4 for the acceptable range of Storage and Transportation environmental conditions.

Always ensure that the unit and accessories are completely dry and free of any condensation before being packed.

## Certificate of Decontamination

We respect the health and safety of our clients and employees, and request that any products or accessories being returned are decontaminated in accordance with the procedure below.

### 1. Decontamination Procedure

Thoroughly clean all outside surfaces of the product (including any accessories, power cords, manuals, packaging, etc) with a damp cloth soaked with suitable disinfectant solution (such as Virkon).

Allow to dry fully before packing.

### 2. Decontamination Declaration

Company Name: \_\_\_\_\_

Address: \_\_\_\_\_

Product Code: **xxsealer - 953-A VTS**

Serial Number: \_\_\_\_\_

Reason For Return: \_\_\_\_\_

Where Product Used: \_\_\_\_\_

Please tick the appropriate option(s) below:

- I certify that I have decontaminated the product as per the above procedure.

Decontaminant Used: \_\_\_\_\_

- I certify that the product has not been exposed to any chemical or biological materials.

Title: \_\_\_\_\_

Name: \_\_\_\_\_

Signature: \_\_\_\_\_

Date: \_\_\_\_\_

Telephone: \_\_\_\_\_

Email: \_\_\_\_\_



## 9 Warranty and Returns

BJS Biotechnologies Ltd. warrants the **xxsealer** product, when purchased new and installed and operated in accordance with the instructions of this manual, to be free from defects in materials and workmanship, and will repair or replace, at their discretion, any unit or accessory which exhibits such defects.

In no event will BJS Biotechnologies be liable for any indirect, incidental or consequential damages resulting from any defect or warranty claim.

**NOTE**

Unspecified use or unauthorised modification of any part of the **xxsealer** unit or its accessories or the use or attachment of any adaptor or peripheral not supplied, specified or sanctioned by BJS Biotechnologies will invalidate this warranty.

This warranty is provided to the original purchaser of the product for one year from the date of purchase.

Under the terms of this warranty, the product must be returned in its original packaging, transportation prepaid by the sender, with a copy of the Proof of Purchase and a detailed description of the problem.

**WARNING**

The product must be decontaminated using the procedure detailed in Section 8.4 and a Certificate of Decontamination supplied with any return. If the product is considered too hazardous to be shipped, please contact BJS Biotechnologies on the number given on Page 2 of this manual for further instructions.

Please contact your distributor (or BJS Biotechnologies on the number given on Page 2 of this manual) to assess the defect and arrange the return the product if necessary.

## 10 Product Disposal

At end-of-life, this product must be disposed of in accordance with your local authority regulations for the disposal of potentially hazardous waste and electronic equipment.

The unit and its accessories should be decontaminated using the procedure detailed in Section 8.4 before disposal or shipping.



Do not dispose of this product into unsorted municipal waste or public landfill.

Please contact your distributor (or BJS Biotechnologies at the address on Page 2 of this manual) for details of how to correctly dispose of this product.