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TEMPO® QC*For microbiological control only*

Kit for the verification of TEMPO Reader measurement performance

SUMMARY AND EXPLANATION

TEMPO QC should only be used with the TEMPO system. The TEMPO QC kit is used to check TEMPO Reader measurement performance. Each control kit is intended for single use only.

The kit should be used to confirm that the instrument is functioning correctly and ensure that the fluorescence readings are accurate.

PRINCIPLE

The TEMPO QC kit consists of 5 vials of standard solution and 5 control cards. All the kit components are ready-to-use.

Each vial and each card is identified by a pictogram (8 or 20).

Each vial is associated with a card which has the same pictogram. The TEMPO Filler fills each card with the solution contained in the matching vial. It also seals the card to avoid any risk of contamination during subsequent handling.

The solution contained in the cards emits a calibrated fluorescent signal which depends on its concentration of 4-Methylumbelliferone (8 or 20 $\mu\text{mol.l}^{-1}$). The light signal detected by the TEMPO Reader, is evaluated by comparison with an expected value.

CONTENT OF THE KIT (Kit for 1 control):

- 3 TEMPO QC 13 ml vials.8 μM
- 2 TEMPO QC 13 ml vials.20 μM
- 3 TEMPO QC cards.8 μM
- 2 TEMPO QC cards.20 μM
- 1 package insert

COMPOSITION OF THE VIALS OF STANDARDS

	4-Methylumbelliferone
Standard 8 μM	8 $\mu\text{mol.l}^{-1}$
Standard 20 μM	20 $\mu\text{mol.l}^{-1}$

Volume of the vials: 13 ml

WARNINGS AND PRECAUTIONS

- **For microbiological control only.**
- **For professional use only.**
- This kit is for single use only.
- Do not incubate the control cards.
- It is critical to correctly match the vials and control cards with the same pictograms (8 or 20).
- Do not use kit components after the expiry date indicated on the box label.
- Do not mix components from different lots.
- Before use, check that the packaging and components are intact.
- Allow the components to come to room temperature in their box before use.
- Do not use visibly deteriorated cards.
- The TEMPO Reader, TEMPO Filler and racks should be regularly cleaned and decontaminated (see the User's Manuals).

- Any change or modification in the procedure may affect the results and must be validated by the laboratory. bioMérieux will not be held liable for results obtained following any changes or modifications in procedures not validated by bioMérieux. In addition, such changes or modifications may void all warranties.

STORAGE CONDITIONS

- Store the components of the TEMPO QC kit between 2°C and 8°C in their box protected from the light.
- If stored according to the recommended conditions, all components are stable until the expiry date indicated on their label.
- After opening the kit, do not leave the components exposed to direct light.

TEST FREQUENCY

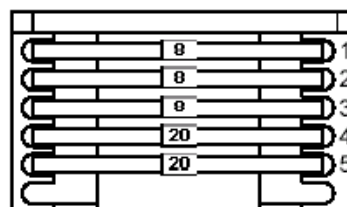
It is recommended to check the TEMPO Reader measurement performance once a month.

INSTRUCTIONS FOR USE

For complete instructions, see the TEMPO system User's Manuals (TEMPO Check chapter).

1. Prior to use, allow the TEMPO QC kit to come to room temperature, then remove the components from their box.
2. Check the expiry date of each of the components.
3. Put each vial in the filling rack and insert the cards with the same pictogram opposite the vial. The transfer tube of each card should be placed inside the matching vial.
4. Place the rack in the TEMPO Filler and start the filling-sealing cycle. The TEMPO Filler simultaneously fills the cards, then cuts and seals the transfer tubes.
5. Remove the filling rack from the TEMPO Filler.

NOTE: There is a residual volume of standard solution in the vials after the filling cycle.
6. Take the cards out of the filling rack and transfer them into the reading rack, ensuring that the label of each card is facing the user (towards the rack handle) and that the cards are properly inserted into the following positions:
 - the three 8 μM cards: rack positions 1, 2 and 3.
 - the two 20 μM cards: rack positions 4 and 5.

7. Dispose of the used vials and transfer tubes into an appropriate receptacle.

READING AND INTERPRETATION

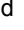
1. Open the TEMPO Check application on the reading station.
2. Introduce the reading rack into the TEMPO Reader and start the reading. The TEMPO Reader checks that the cards are in the correct positions in the rack, by scanning each of the barcodes, then detects and interprets the fluorescence for each card.
3. Consult the test results display and management screen. Once the reading is completed, the results are automatically analyzed by the computer, which determines whether the reader is functioning correctly, by comparing the results to the range of expected fluorescence values.
4. Remove the cards from the rack and dispose of them into an appropriate receptacle.

EXPECTED RESULTS

When reading of the TEMPO QC cards is completed, the screen for the display and management of the results of the TEMPO Reader measurement performance controls displays the results:

- : The reader functions correctly.
- : The reader functions correctly but you must be extremely vigilant when the next controls are performed.
- : See "Troubleshooting" section below.

Troubleshooting

- Check if the storage conditions (2-8°C and no exposure to direct light) and instructions for use (allow the kit to come to room temperature before use) have been respected.
- Change the position of the card with  status * and read again.
- If the result of the control is still red () , perform the control again using another QC kit which has been correctly stored and allowed to come to room temperature before use.
- If the result of the control obtained with this second QC kit is still red () , change the position of the card with  status * and read again.
- If the result of the control is still red () , please contact bioMérieux SA or its local representative.

* making sure the 8µM cards are in positions 1, 2 and 3 and the 20 µM cards are in positions 4 and 5.

NOTE: A maximum number of 2 readings can be performed with each kit.

WASTE DISPOSAL

It is the responsibility of each laboratory to handle waste and effluents produced according to their nature and degree of hazardousness and to treat and dispose of them (or have them treated and disposed of) in accordance with any applicable regulations.

INDEX OF SYMBOLS

Symbol	Meaning
	GB: Catalogue number US: Catalog number
	Manufacturer
	Temperature limitation
	Protect from light
	Use by
	Batch code
	Consult Instructions for Use

WARRANTY

bioMérieux disclaims all warranties, express or implied, including any implied warranties of MERCHANTABILITY AND FITNESS FOR A PARTICULAR USE. bioMérieux shall not be liable for any incidental or consequential damages. IN NO EVENT SHALL BIOMERIEUX'S LIABILITY TO CUSTOMER UNDER ANY CLAIM EXCEED A REFUND OF THE AMOUNT PAID TO BIOMERIEUX FOR THE PRODUCT OR SERVICE WHICH IS THE SUBJECT OF THE CLAIM.

 **bioMérieux® SA**
 au capital de 12 029 370 €
 673 620 399 RCS LYON
 69280 Marcy-l'Etoile / France
 Tel. 33 (0)4 78 87 20 00
 Fax 33 (0)4 78 87 20 90
<http://www.biomerieux.com>

bioMérieux, Inc
 Box 15969,
 Durham, NC 27704-0969 / USA
 Tel. (1) 919 620 20 00
 Fax (1) 919 620 22 11

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Brasil: Distribuído por bioMérieux Brasil, S.A. - Estrada do Mapuá, 491 - Jacarepaguá - R.J. - CEP 22710-261