IsoTherm-System®

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eppendorf

Copyright© 2008 Eppendorf AG, Hamburg. No part of this publication may be reproduced without the prior permission of the copyright owner.

Trademarks

eppendorf® and eppendorf Thermomixer® are registered trademarks of Eppendorf AG, Hamburg, Germany.

Thermomixer™ is a trademark of Eppendorf AG, Hamburg, Germany.

Trademarks are not marked in all cases with ™ or ® in this manual.

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
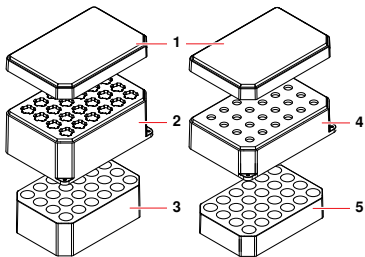
1.1 Main illustration

IsoRack and IsoPack

1 Tube holder cover	2 IsoRack 1.5/2.0 ml for 24 1.5/2.0 ml Safe-Lock tubes
3 IsoPack 1.5/2.0 ml for 1.5/2.0 ml IsoRack	4 IsoRack 0.5 ml for 24 0.5ml Safe-Lock tubes
5 IsoPack 0.5ml for 0.5ml IsoRack	

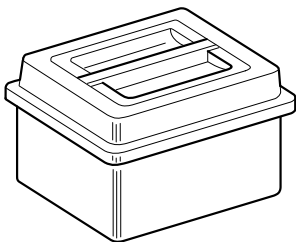IsoSafe

1.2 Features

With the IsoTherm system you can cool samples efficiently and at a constant temperature for several hours and then thaw them gently.

The tube holder further serves as an aid when filling micro test tubes and for provisioning, sorting, distributing, transporting, storing and freezing samples in micro test tubes.

The IsoTherm system includes the components IsoRack, IsoPack and IsoSafe.

Component	Variants	Features
IsoRack (tube holder)	<ul style="list-style-type: none"> • for 0.5 ml Safe-Lock tubes, or • for 1.5/2.0 ml Safe-Lock tubes, each with cover. 	<ul style="list-style-type: none"> • base of an MTP • accepts 24 micro test tube; numbered per row • temperature resistant between -200 and $+100$ °C • UV resistant and suitable for autoclaving • suitable for centrifuging in the MTP rotor with IsoRack adapter ¹ • stackable with or without cover • bore holes at the distances of the MTP grid, allowing for micro test tubes to be filled using multi-channel pipettes • cover fits twist-proof, locks to the tube holder • cover can also be used as base for cool pack and tube holder • floatable with cover (thawing in a water bath)
IsoPack (cool pack)	<ul style="list-style-type: none"> • for 0.5-ml IsoRack or • for 1.5/2.0-ml IsoRack 	both sizes are available in two versions each: <ul style="list-style-type: none"> • IsoPack white for 0 °C • IsoPack blue for -21 °C
IsoSafe (insulating box)	—	<ul style="list-style-type: none"> • polystyrene box with cover for one IsoPack each with corresponding IsoRack • ensures the cooling of the samples

¹ Generally the centrifuging stability depends on the centrifuge, rotor, rotor buckets, adapter, maximum rpm / maximum field of gravity (rcf), operating temperature, centrifugation times, density of the centrifuged liquid, organic solvents, accumulated total weight and proper operation.

1.2.1 Dwell time

For the cool packs 0.5 ml or 1.5/2.0 ml the dwell time (IsoTherm time) has been calculated for 3 use cases each (see measurement in the table):


- Insulating box with cover (a)
- Open insulating box (b)
- Cool pack and tube holder on the tube holder cover (c)

Use case ¹	a		b		c	
	0 °C to 4 °C (white)	-21 °C to 18 °C (blue)	0 °C to 4 °C (white)	-21 °C to -18 °C (blue)	0 °C to 4 °C (white)	-21 °C to -18 °C (blue)
IsoTherm time	>6 h	>3 h	>4 h	>1 to 2 h	> 1.5 to 3 h	> 0.5 to 1 h ²

¹ Framework conditions for measuring the IsoTherm time:

- Maximum filling of the tube holder.
- Micro test tubes filled to the maximum.
- Environment at room temperature, system at rest.
- Measurement of the sample temperature.

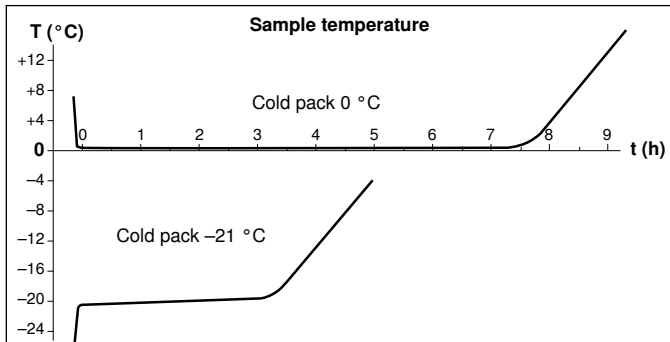² For IsoPack 1.5/2.0 ml: only valid for central bore holes.

2 Intended use

The IsoTherm System is intended exclusively for in-door use and for the cooling of tubes. Use only Eppendorf accessories or accessories recommended by Eppendorf. The IsoTherm system must only be operated by appropriately qualified and instructed personnel. Carefully read the operating manual before working with the IsoTherm system and familiarize yourself with its use.

3 Operation

Caution! Damage of the IsoPack from sharp edges.

Sharp edged components can damage the film on the underside of the IsoPack.

- ▶ Do not use sharp edged components near the IsoPack.
-

Recommendations for pre-cooling

- Cool pack 0 °C: approx. 2 hours (** freezer) or over night (* freezer)
- Cool pack -21 °C: approx. 4 hours at -35 °C (*** freezer, setting: rapid freeze)

Operation

1. Insert pre-cooled cool box into the insulating box.
2. Place tube holder with samples onto the cool box.
3. Close insulating box with a cover.

This prevents the formation of frost.

The samples can be temperature controlled for several hours at 0 to -21 °C.

4 Maintenance

4.1 Cleaning

Clean the IsoTherm system components immediately if sample fluid enters the bore holes or the surface.

1. Clean the components of the IsoTherm system with a mild soap solution.
2. Rinse thoroughly with distilled water.

The IsoRack is UV resistant and can be autoclaved at 121 °C for 20 minutes. The tube holder and cover are resistant to alcohol.

Do not use organic solvents such as phenol, chloroform and acetone.

4.2 Disinfection / decontamination

Warning! Health hazard from contaminated accessories.

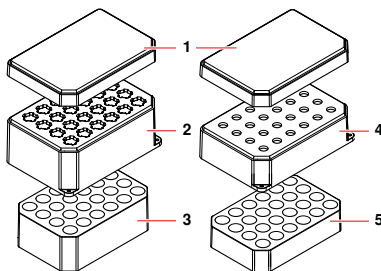
1. Follow the instructions in the decontamination certificate. It is available in PDF format on our homepage (www.eppendorf.com/decontamination).
 2. Decontaminate all the parts you want to dispatch.
 3. Include the fully completed decontamination certificate for returned goods in the package.
-

- ▶ Select a disinfection method complying with the statutory rules and regulations for your area of application. Use e.g. alcohol (ethanol, isopropanol) or alcohol-containing disinfectants.

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Eppendorf AG · 22331 Hamburg · Germany · Tel: +49 40 538 01-0 · Fax: +49 40 538 01-556
Eppendorf North America, Inc. · 102 Motor Parkway, Suite 410 · Hauppauge, NY 11788-5178 · USA