

RNA and DNA purification products from MACHERY-NAGEL

RNA and DNA isolation for veterinary diagnostics

One kit for all isolations

NucleoMag[®] VET



Purification of viral RNA/DNA and bacterial DNA

Magnetic-bead based isolation

Highly sensitive and reliable downstream applications

... one kit for common veterinary sample materials

MACHERY-NAGEL

www.mn-net.com



Since 1911

NucleoMag[®] VET

NucleoMag[®] VET

For purification of viral nucleic acids and bacterial DNA

Select sample-specific pre-lysis procedure

Nucleic acid isolation with universal NucleoMag[®] VET protocol


Downstream detection system

The NucleoMag[®] VET kit provides purification of viral nucleic acids and bacterial DNA from common veterinary samples. Nucleic acids are easily purified from samples such as whole blood, serum, plasma, tissue homogenates, ear notches, and swabs. A sample-specific pre-lysis procedure is recommended followed by an universal purification protocol based on magnetic-bead technology.

Product at a glance

Technology	Magnetic-bead technology
Processing	Automated or manual
Sample material	Animal whole blood and plasma/serum (200 µL), ear notches, tissue (10–30 mg), swabs, feces (500 µL)
Fragment size	300 bp–approx. 50 kbp
Elution volume	50–100 µL

Application data

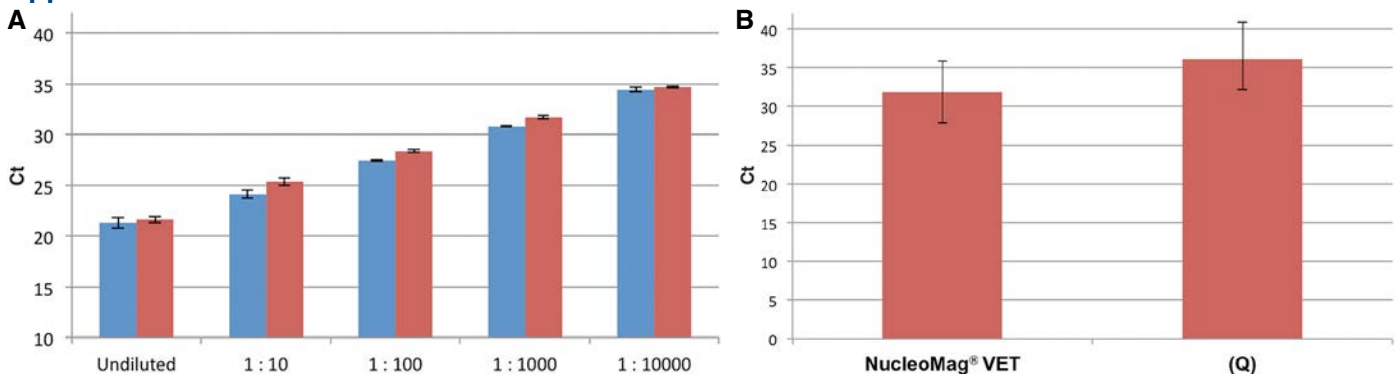

Sensitive detection of viral RNA and DNA from animal samples.

(A) For purifications, serum samples from PRRSV (blue) and PCV2 (red) positive animal samples were diluted 10-fold. Viral RNA (PRRSV) or viral DNA (PCV2) were isolated from 200 µL blood serum sample using NucleoMag[®] VET kit. Purified RNA (PRRSV) was analyzed using Real-time Multiplex RT-PCR (VIROTYPE PRRSV, LDL). Purified DNA (PCV2) was analyzed using Real-time PCR (PCV2 Primer). *Data was generated in co-operation IVD GmbH Hanover, Germany.*

(B) PCV2 positive animal tissue samples (n=22, 20 mg each) were used for viral DNA purification using indicated kits. Purified viral DNA was analyzed by Real-Time PCR (PCV2 Primer). Out of 22 samples 20 (MN) or 14 (Q) samples were tested positive with Ct mean values shown in the graph. *Data kindly provided by Italian customer.*

Ordering information

Product	Preps	REF
NucleoMag [®] VET	1 x 96 / 4 x 96	744200.1 / .4

Your local distributor

www.mn-net.com

MACHEREY-NAGEL



MACHEREY-NAGEL GmbH & Co. KG · Neumann-Neander-Str. 6–8 · 52355 Düren · Germany

Germany and international:
Tel.: +49 24 21 969-0
Fax: +49 24 21 969-199
E-mail: info@mn-net.com

Switzerland:
MACHEREY-NAGEL AG
Tel.: +41 62 388 55 00
Fax: +41 62 388 55 05
E-mail: sales-ch@mn-net.com

France:
MACHEREY-NAGEL EURL
Tel.: +33 388 68 22 68
Fax: +33 388 51 76 88
E-mail: sales-fr@mn-net.com

USA:
MACHEREY-NAGEL Inc.
Tel.: +1 484 821 0984
Fax: +1 484 821 1272
E-mail: sales-us@mn-net.com

