

Total Protein Quantification from Placenta using the Bead Ruptor Elite™ Bead Mill Homogenizer for Sample Preparation

Gabriella Ryan, M.S.; Rodney J Nash, Ph.D.
OMNI International, a PerkinElmer company

Summary

The impact of proteomics research on placenta tissue has allowed scientists to discover patterns in key biomarkers related to both placental and gestational syndromes and pathologies¹. Specifically, at the bench, uncovering changes in biomarkers related to preeclampsia², gestational diabetes³, and TORCH⁴ (Toxoplasmosis, "Others", Rubella, Cytomegalovirus, Herpes simplex virus) syndrome, to name a few, are propelling researchers and clinicians forward in being able to diagnose and treat these conditions during pregnancy.

Upstream of protein-based assays like Western blot, ELISA, LC-MS/MS is a sample preparation step involving homogenization of fixed placenta tissue, which lyses cells and releases analytes of interest. Standard homogenization methods involving harsh detergents and enzymes can be time consuming and ineffective procedures for tough samples like placenta. Bead mill homogenization on the Bead Ruptor Elite™ Bead Mill Homogenizer is a quick and effective alternative to standard chemical-based sample preparation methodologies, while still producing a homogenate suitable for downstream analysis.

Herein, we outline sample preparation of fixed placenta tissue on the Bead Ruptor Elite for total protein quantification.

Materials and Methods

Equipment

- Bead Ruptor Elite™ Bead Mill Homogenizer (Cat # 19-042E)
- Bead Ruptor Elite™ Bead Mill Homogenizer 2 mL Tube Carriage (Cat # 19-373)
- Bead Ruptor Elite™ Bead Mill Homogenizer Finger Plate (Cat # 19-370)
- 2 mL Hard Tissue Homogenizing Mix (Cat # 19-628)



**2 mL Hard Tissue
Homogenizing
Mix**



**Bead Ruptor Elite™
Bead Mill Homogenizer**



**Bead Ruptor Elite™ Bead
Mill Homogenizer Finger Plate**



2 mL Tube Carriage Kit



OMNI
International
a PerkinElmer company

Procedure

Sample Acquisition

De-identified paraformaldehyde-fixed human placenta tissue samples were obtained from Emory University Hospital, Tissue Procurement Lab under standard IRB protocols.

Total Protein BCA Assay

200 mg of placenta tissue was weighed out and transferred to a 2 mL Hard Tissue Homogenizing Mix Tube (Omni International, Cat No. 19-628) along with 600 μ L of phosphate buffered saline (PBS) (Gibco, Cat No. 20012027). Samples were weighed out with a tolerance \pm 10 mg. Placenta samples were processed on the Bead Ruptor Elite™ Bead Mill Homogenizer at 5 m/s for 2 cycles of 30 seconds (Table 1). Homogenate was then transferred to a microplate and prepared according to the instructions for 'Microplate Procedure' in the BCA Protein Assay Kit (Thermo Scientific, Cat No. 23225). Working Reagent and Bovine Serum Albumin (BSA) standards A-G (Table 2) were also prepared according to dilution scheme provided in the kit instructions. Prepared microplate sample absorbance was measured at 562 nm, along with included BSA Standards. Blank 562 nm absorbances were used for further construction of standard curve and analysis of measured protein concentration.

Sample Type	Sample Weight (mg)	Speed (m/s)	Time (sec)	Cycles	Dwell Time (sec)
Placenta	200	5	30	2	10

Table 1: Placenta Sample Homogenization Summary.

Results

A complete homogenate was obtained after 1 minute of processing time (Figure 1). Downstream BCA Total Protein quantification yielded an average protein concentration of 1064 μ g/mL (Table 3). Protein concentration was determined based on the equation generated from the standard curve, using BSA standards for measurement of absorbance at 562 nm.



Figure 1: Pre and post-processing images of 200 mg placenta tissue.

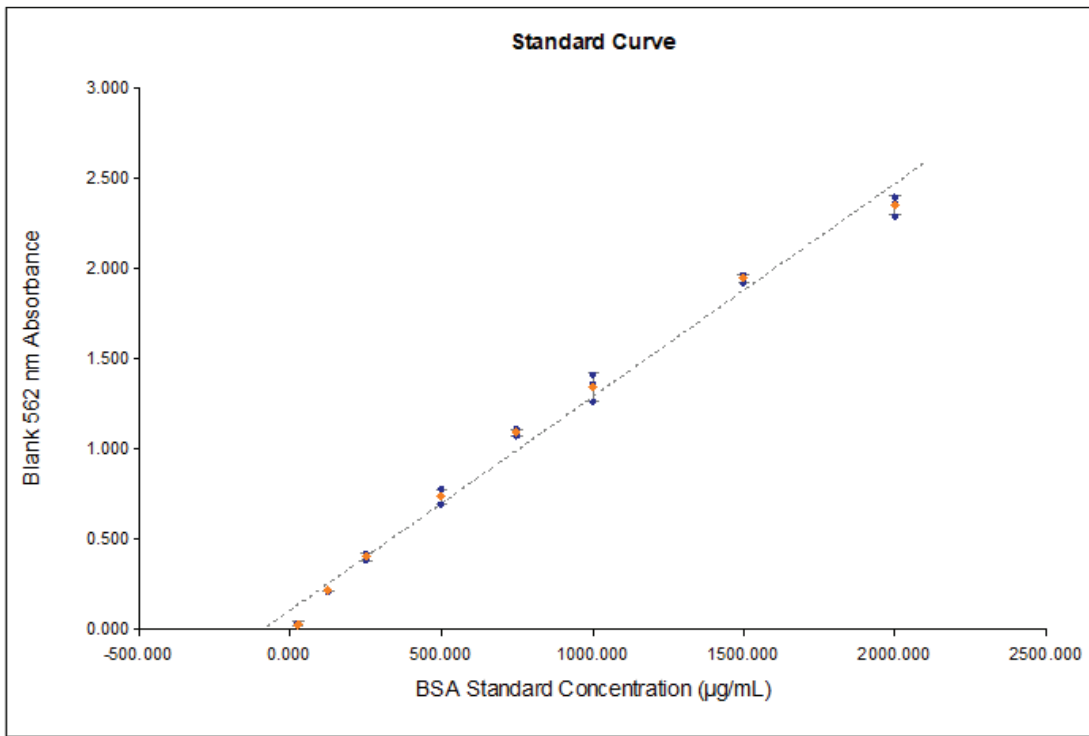


Figure 2. Standard curve generated using BSA standards A-G from the BCA Protein Assay Kit. Equation generated from standard curve: $y=mx + b$, where y is defined as the Blank 562 nm measurement, m is defined as 0.00118, b is defined as 0.104 and x is protein concentration ($\mu\text{g/mL}$).

Standard	Known Concentration ($\mu\text{g/mL}$)	Average Measured Concentration ($\mu\text{g/mL}$)
A	2000	1904
B	1500	1562
C	1000	1051
D	750	837
E	500	536
F	250	235
G	125	92
PBS	0	0

Table 2. Average BSA Standard Protein Concentration between prepared triplicate standards.

Sample	Calculated Protein Concentration (µg/mL)
1	938
2	1129
3	1123

Table 3. Calculated protein concentrations from homogenized placenta tissue.

Conclusions

The Bead Ruptor Elite™ Bead Mill Homogenizer along with 2 mL Hard Tissue Homogenizing Mix offers scientists a sample preparation solution for placenta tissues yielding a complete homogenate compatible for downstream total protein quantification. As seen in the above experiments, high yield protein concentration recovered from homogenized placenta is a suitability indicator for further proteomics assays like ELISA, LC-MS/MS and Western blot.

References (AMA)

1. Robinson JM, Ackerman WE 4th, Kniss DA, Takizawa T, Vandr  DD. Proteomics of the human placenta: promises and realities. *Placenta*. 2008;29(2):135-143. doi:10.1016/j.placenta.2007.12.005
2. Chang X, Bian Y, Wu Y, Huang Y, Wang K, Duan T. Endocan of the maternal placenta tissue is increased in pre-eclampsia. *Int J Clin Exp Pathol*. 2015;8(11):14733-14740. Published 2015 Nov 1.
3. Burlina S, Banfi C, Brioschi M, et al. Is the placental proteome impaired in well-controlled gestational diabetes?. *J Mass Spectrom*. 2019;54(4):359-365. doi:10.1002/jms.4336
4. Macedo-da-Silva J, Marinho CRF, Palmisano G, Rosa-Fernandes L. Lights and Shadows of TORCH Infection Proteomics. *Genes (Basel)*. 2020;11(8):894. Published 2020 Aug 5. doi:10.3390/genes11080894

**For research use only.
Not for use in diagnostic procedures.**

Ordering Information

Contact Your Local OMNI International Representative for:

- Bead Ruptor Elite™ Bead Mill Homogenizer (Cat # 19-042E)
- Bead Ruptor Elite™ Bead Mill Homogenizer 2 mL Tube Carriage (Cat # 19-373)
- Bead Ruptor Elite™ Bead Mill Homogenizer Finger Plate (Ca t# 19-370)
- 2 mL Hard Tissue Homogenizing Mix (Cat # 19-628)



Contact Info for your Local OMNI International Representative:

sales@OMNI-inc.com

or

<https://us.omni-inc.com/contact-us>



OMNI International, a PerkinElmer company
935 Cobb Place Blvd NW
Kennesaw, GA 30144
United States
770.421.0058 | 800.766.4431